

caArray 022 - Changing the ownership of the experiment in caArray

Question: Can I change the ownership of the experiment in caArray?

Topic: caArray Usage

Release: caArray 2.2.1 and up

Date entered: 06/04/2009

Details about the Question

I have created a new experiment for the investigator. Now I need to transfer the ownership of the experiment to my investigator. Can I do that?

Answer

Yes. Changing ownership is a new feature implemented in caArray Version 2.2.1 (Released May, 2009). The ownership transfer can only be performed by a caArray user with system administrator privileges.

Upon login as system administrator, a link called "Manage Ownership" will appear under "User management" in the left menu. To Change the ownership, please follow steps below:

1. System administrator (SA) clicks **"Manage Ownership"**.
2. After providing First Name or Last Name, SA clicks "Filter". A list of users who meet the filtering criteria is returned.
3. SA selects the Username of the user who currently owns the experiment, and clicks "Find Assets" to find the experiments associated with this user.

Finding the owner and the owner's assets

WELCOME TO CAARRAY

HOME

BROWSE

EXPERIMENT MANAGEMENT

MY EXPERIMENT WORKSPACE

CREATE/PROPOSE EXPERIMENT

USER MANAGEMENT

MANAGE USERS

MANAGE OWNERSHIP

MANAGE COLLABORATION GROUPS

CURATION

MANAGE ARRAY DESIGNS

MANAGE PROTOCOLS

MANAGE VOCABULARY

ABOUT CAARRAY

WHAT IS CAARRAY?

RELEASE NOTES

TECHNICAL DOCUMENTATION

TRAINING

HELP

GLOBAL QUICK LINKS

NATIONAL CANCER INSTITUTE (NCI)

NCI CENTER FOR BIOINFORMATICS (NCIB)

Manage Ownership

Users

Search for users by choosing filter criteria. Select the user who's assets you want to re-assign, then press the Find Assets ... button.

1

Last Name:

First Name:

Organization:

Clear Filter

14 items found, displaying all items.

	Username	Member Name	Organ
☐	caarraytest	Tester_caArray	
☐	bjagla	Yagla, Bernd	CUMC
☐	zli	Li, Zhong	CUMC
☒	flin	Lin, Fan	CUMC
☐	cu_reviewers	read-only reviewer	
☐	ksmith	Smith, Ken	
☐	caarrayadmin	Administrator_caArray	
☐	caarrayuser	User_caArray	

Find Assets ... 2

4. SA selects the experiment that needs a new owner.

- SA clicks on "Find New Owner".
- Select experiment and finding the new owner

Assets owned by Fan Lin

Select the assets to re-assign, then press the "Find New Owner..." button below to select the new owner.

Experiments	Experiment ID	Experiment Title	Assay Type
☐	lin-00010	Transcriptional profiling of a...	Gene Expression
☐	lin-00011	tl_test02	Gene Expression
☐	lin-00014	Example Multi SDRF Files	Gene Expression
☐	lin-00015	manual enter annotation data	Gene Expression
☐	lin-00018	Extra file test	Gene Expression
☐	lin-00020	test1	Gene Expression
☐	lin-00025	test	Gene Expression
☐	lin-00026	SNP test	SNP
☒	lin-00031	Epidemioid Carcinoma	Gene Expression

Collaboration Groups

2 items found, displaying all items.

Collaboration Group Name	Group Members
caarraytest	caArray User, caArray Tester
Reviewer	reviewer read-only

Find New Owner...

- SA selects the New Owner
 - SA clicks "Assign"
- The owner of this experiment is now changed to the new owner as shown in the image.

Assign New Owner to the experiment

Last name:

First name:

Organization:

13 items found, displaying all items.

Username	Member Name
☒ caarraytest	Tester, caArray
☐ bjagla	Yagla, Bernd
☐ zli	Li, Zhong
☐ cu_reviewers	read-only, reviewer
☐ ksmith	Smith, Ken
☐ caarrayadmin	Administrator, caArray
☐ caarrayuser	User, caArray
☐ researchscientist	ResearchScientist, ResearchScientist
☐ labadministrator	LabAdministrator, LabAdministrator
☐ labscientist	LabScientist, LabScientist
☐ biostatistician	Biostatistician, Biostatistician
☐ systemadministrator	SystemAdministrator, SystemAdministrator
☐ collaborator	Collaborator, Collaborator

Assign

Have a comment?

Please leave your comment in the [caArray End User Forum](#).